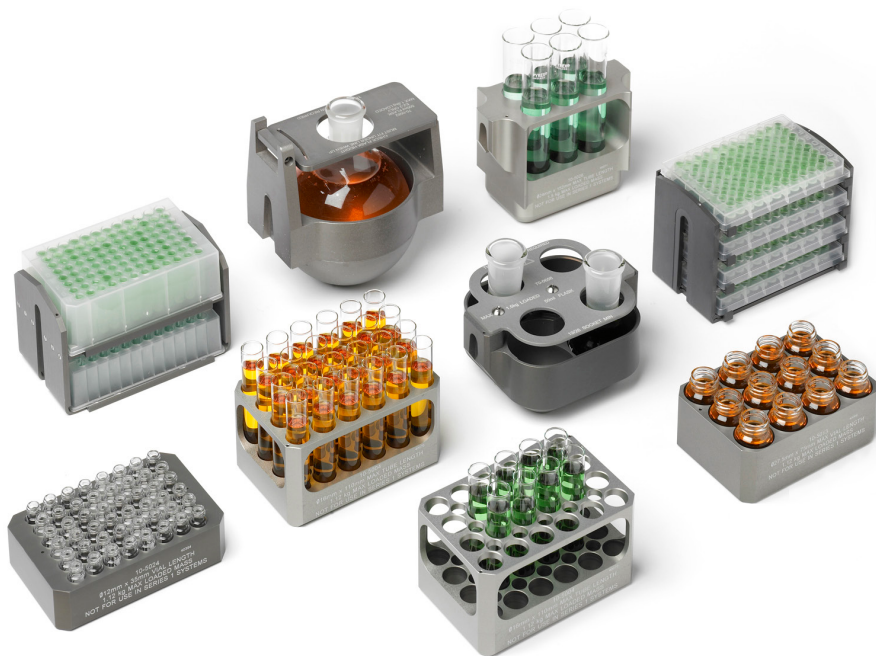




Accessories

for Series 3 & EZ-2 Evaporators



Part Number 10-1478 - Revision 8.5 - February 2022

Disclaimer

Copyright © 2021 Genevac Ltd.

All trademarks referenced in this document are registered or common law trademarks of SP Industries, Inc. or its affiliates in the U.S. and/or other countries.

Original Instructions

The U.S. English version of this document is the original instructions.

All other languages are a translation of the original instructions.

Genevac Ltd.

The Sovereign Centre

Farthing Road

Ipswich, Suffolk

IP1 5AP, United Kingdom

+44 (0) 1473 240000

SP Service (US) +1 (877) 548-4666

SP Service (OUS) +44 (0)1473 243000

SP Service Email: service@spscientific.com

Website: scientificproducts.com

This Accessories Manual contains confidential and proprietary information of Genevac Ltd. (hereinafter “Company”) and may be used only by a recipient designated by and for purposes specified by Company. You may not disclose the contents of this Accessories Manual to any unauthorized third party. You agree that you will only provide this Accessories Manual to officers, employees, contractors, or other service providers who have a need to know this information and who are bound by written confidentiality obligations.

You may only use this copy of this Accessories Manual. Unless otherwise authorized by Company in writing, you have no right to reproduce, distribute, modify, create derivative works, publicly display or perform this Accessories Manual. Permitted copies of this document must retain all proprietary notices contained in the original.

Your purchase of the equipment is governed by Company’s then-current terms of purchase. The information in this document is subject to change at any time in Company’s sole discretion without notice. As such, we suggest that you regularly confirm with Company that you are using the most current version of this document. Company is free to modify any of its products and services, in any manner and at any time.

EXCEPT AS EXPRESSLY PROVIDED HEREIN, THE CONTENTS OF THIS DOCUMENT SHALL NOT CONSTITUTE ANY WARRANTY OF ANY KIND, EITHER EXPRESSED OR IMPLIED, INCLUDING BUT NOT LIMITED TO THE IMPLIED WARRANTIES OF MERCHANTABILITY AND/OR FITNESS FOR A PARTICULAR PURPOSE OR GIVE RISE TO ANY LIABILITY OF COMPANY, ITS AFFILIATES OR ITS SUPPLIERS.

Table of Contents

Introduction	4	Bottle Holders	28
Nomenclature	4	Flask Holders	29
Sample Holder Compatibility	4	SampleGenie™	31
Sample Holder Types	5	60 ml SampleGenie	32
Heat Transfer	6	60 ml Drying Adapters	33
Heat-transfer Plates for Microplates	6	60 ml Lyophilisation/Concentration Adapters	33
Thermal Cushions & Tip Cushions	7	125 ml SampleGenie	34
Protective Coatings	7	125 ml Drying Adapters	35
Sample Holders for Special Applications	7	125 ml Lyophilisation/Concentration Adapters	35
Minimum Order Quantities		250 ml SampleGenie	36
First-time Use	8	250 ml Drying Adapters	37
Maintenance	8	250 ml Lyophilisation/Concentration Adapters	37
Inspection	8	Miscellaneous	38
Cleaning	8	Infinity Trolley	38
Glassware	9	Heat-Transfer Plates for Microplates	40
Labelling of Glassware	9	Gilson and Waters Fraction Collection Racks	41
Swings	10	Thermal Cushions	43
Side-bridge Swing	10	Tip Cushions	45
Fast-Stack Swing - Shallow Well	10	Viton Cushion	45
Fast-Stack Swing - Deep Well	10	Spare Parts	46
Sample Holders	11	EZ-2 Consumables	46
Tube holders for Side-bridge Swings	11	EZ-2 Mk2 & Mk3	46
One-Piece Tube Holders	16	EZ-2 4.0	47
Vial-Holders for Side-Bridge Swings	19	Series 3 Consumables	49
One-Piece Vial Holders	25	Series 3 SampleGuard Probes	51
Centrifuge Tube Holders	26	Genevac Patents	55
Microcentrifuge Tube Holders	27		

Introduction

SP Genevac sample holders are manufactured from high-grade, solid aluminium to exacting specifications, including precise mass-balancing to ensure smooth operation of the evaporator.

Selection of the correct sample holder is a critical factor in optimising evaporation performance for your selected application. If you have any questions regarding the best sample holder for your application, please contact your local SP representative.

The common holders and SampleGenie adaptors are listed here. For the full range, please contact you local SP representative.

Nomenclature

Throughout this brochure, the following terms are used:

Tube Holder: A sample holder that accommodates tubes with rounded ends.

Vial Holder: A sample holder that accommodates vials with flat ends.

Bottle Holder: A sample holder that accommodates bottles with flat ends.

Flask Holder: A sample holder that accommodates round-bottomed flasks.

Sample holders intended for use with tubes with conical or tapered ends are identified individually.

Sample Holder Compatibility

Sample holders within this brochure are intended for use in SP Genevac Series 3, Series 2 and EZ-2 series centrifugal evaporators. They are not suitable for use in other evaporators or similar products.

Sample holder compatibility is indicated as follows:

Compatibility	
EZ-2	✓
S2	✓
S3	✓

Some sample holders are subject to limitations of use, or require accessories to function. Limitations and/or required accessories will be indicated as follows:

NOTE: Requires Cushion (supplied)

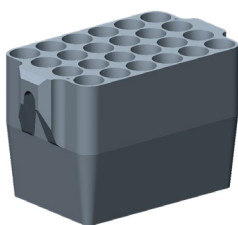
Sample Holder Types

A comprehensive range of sample holders is available to suit tubes, vials, beakers, flasks, microplates etc. Sample holders come in two varieties: one-piece holders, and two-part holders with side-bridge swings.

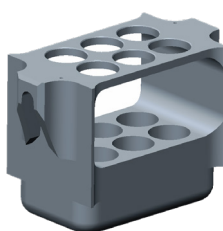
NOTE: The maximum weight of the swing, holder and sample shall not exceed 1.5Kg.

One-piece Sample Holders

One-piece sample holders are those in which the swing is integral to the holder.



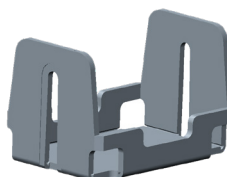
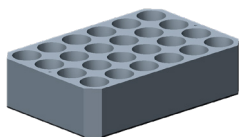
One-piece sample holder



*One-piece sample holder with
integral top-plate*

Two-part Sample Holders

Two-part sample holders are those that consist of a sample holder and separate side-bridge swing.



Separate sample holder and side-bridge swing

All sample-holders are designed to lift cleanly out of the rotor to allow for easy sample loading.

Heat Transfer

A major factor in determining the speed of evaporation is the rate at which heat is transferred into the sample. Greater heat transfer means faster evaporation. In SP Genevac Series 3 and EZ-2 Series evaporators, heat transfer is optimised via the following methods:

High-quality Aluminium Rotors and Sample Holders

SP Genevac evaporators use high-quality aluminium, ensuring the highest thermal conductivity. This allows the most efficient heat transfer, at the lowest weight.

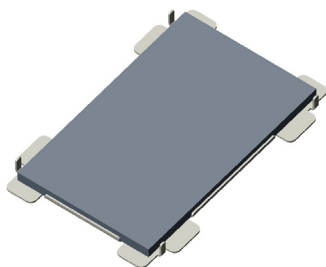
Precision Engineering

Sample holder tube wells must be manufactured to high tolerance to ensure no gaps between the wall and the glassware that would reduce heat transfer or result in breakage. Centre of gravity and mass must be correct to ensure even heat distribution and rotor balance. Precision engineering combines with over two decades of experience to ensure perfect evaporation every time.

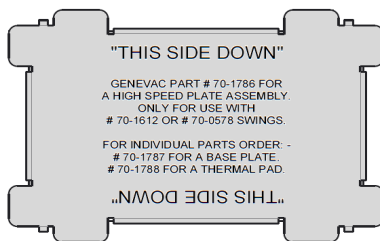
Heat-transfer Plates for Microplates

Heat-transfer plates provide improved evaporation results when using microplates. Improvements will be seen in evenness of sample drying across the plate, and for some applications, heat-transfer plates may help to reduce evaporation times by up to 50%. Improvements are most notable when using deep-well or mid-well plates.

When using heat-transfer plates with pliable thermal pads, the plate should be inserted into the swing with the metal base-plate facing downwards (in contact with the swing). The microplate is loaded onto the top of the pad.



Heat-transfer plate



Heat-transfer pad base-plate

NOTE: Heat-transfer plates are for use with microplates only.

NOTE: Heat-transfer plates **shall** not be fitted with out microplates.

NOTE: Thermal pads are consumables and be replaced when cracks are observed.

Thermal Cushions & Tip Cushions

Thermal Cushions and Tip Cushions provide improved evaporation results in a similar manner to thermal pads used with heat-transfer plates, and prolong glassware lifetimes by helping to absorb the high centrifugal load applied by SP Genevac evaporators. Use of thermal or tip cushions is mandatory with some sample holders.

Sample holders that require the use of thermal or tip cushions are supplied with a starter pack of cushions to allow immediate use.

Cushions are consumable items and must be replaced when worn or damaged. SP recommends purchasing additional cushions to allow for periodic replacement.

Protective Coatings

Standard aluminium sample holders are hard anodised (water sealed) for protection against all common organic solvents, including strong acids such as TFA, HCl and Thionyl chloride.

NOTE: HCl and Thionyl chloride should only be evaporated in systems that are HCl-compatible. These systems are specifically built to resist attack from HCl and Thionyl chloride.

NOTE: The colouring of the anodised surface finish may vary from batch to batch; this is normal and does not affect the performance of the sample holder.

Sample Holders for Special Applications

As the market leader in centrifugal evaporation, SP manufacture sample holders to suit almost every common application. If you have a special application or unusual tube size that is not available in the current range of sample holders, SP offers a bespoke sample holder design service. Please contact your local SP representative for more information.

Minimum Order Quantities

A minimum order quantity of two applies to sample holders, side-bridge swings, heat-transfer plates and SampleGenie assemblies.

First-time Use

When using sample holders for the first time, always check that they are compatible with the evaporator. Load the sample holders with empty tubes (or vials, plates etc. as appropriate) and place the holder and swing assembly onto the rotor. Swing the rotor by hand, allowing the sample holder to reach its fully horizontal state, to make sure that no part of the assembly touches the rotor or chamber.

Maintenance

Inspect and clean the sample holders on a regular basis, and always following any incident of solvent spillage or glass breakage.

Inspection

Visually inspect the sample holders and swings to check for any spilled solvent or debris, especially in the sample holder wells. Spilled solvent can cause debris to adhere to the sample holders, creating raised points that place additional stress on the glassware.

Superficial surface damage such as scratches do not affect the performance of the sample holder. If deep pits or crevices begin to form, or if there is any structural damage (e.g. bending, deformation), immediately stop using it. Contact the SP Service department for evaluation.

Repeated use of sample holders with strong acids may cause colour fading over time. This is normal and does not affect the integrity of the sample holders.

Check the state of all thermal or tip cushions. Replace cushions at the first sign of wear or damage to reduce the risk of glass breakage.

Cleaning

Clean off any adhered dirt or debris by washing the sample holders in soapy water, or with a cleaning solvent such as methanol or acetone. Once washed, rinse the sample holders in clean water. Alternatively, wash in a dishwasher at 90°C. Once washed, inspect the sample holder again to ensure that all adhered dirt, debris and solvent residue is removed.

Always dry the sample holders thoroughly after cleaning. Never place wet sample holders in the evaporator.

NOTE: The most likely cause of glassware breakage is a small fragment of glass from a previous glass breakage which has become stuck in a sample holder well. It is therefore very important that sample holders are cleaned thoroughly after any occurrence of glass breakage.

Glassware

Some SP Genevac evaporators subject samples to accelerations in excess of 500 times gravity. This allows for excellent evaporation, but places a high degree of stress on the glassware. Selection of suitable glassware is essential to avoid glassware breakage and ensure perfect evaporation.

Always use high-quality glassware from recognised manufacturers of international repute. Experience has shown that 'own-brand' catalogue glassware can have highly variable dimensions and manufacturing quality; this kind of glassware is not recommended for use in SP Genevac evaporators due to excessive risk of breakage.

When selecting glassware, ensure the following:

- Reputable scientific glassware manufacturer
- High quality, machine-made borosilicate glass
- Minimum 1.2 mm wall thickness

Re-use of glassware in centrifugal evaporators gradually reduces the strength of the glassware. Replace all glassware regularly to avoid breakage.

Damaged Glassware

Even a small scratch can reduce the strength of glassware by a significant degree. **Do not use glassware that is scratched, chipped, cracked or otherwise damaged. Do not use metal spatulas, wire brushes or similar hard instruments to clean glassware or remove dried samples.**

Labelling of Glassware

There are two common methods of labelling glassware in laboratory environments. Both types can cause problems in the evaporator.

Paper/Thin Film Labels

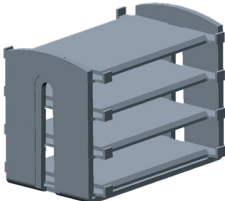
The label effectively increases the diameter of the tube (especially if it wraps around the tube more than once). When the tube is not fully snug in the holder, heat transfer is reduced and any raised points (e.g. tiny fragments of glass from a previous breakage, small lumps of paper towel left behind after cleaning, or the label itself) place additional stress on the glassware.

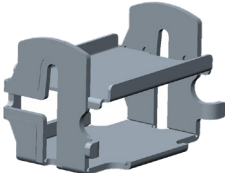
Ceramic Etched

Ceramic-etched labelling weakens the glass. SP recommend testing new tubes for breakage before purchasing glassware in bulk.

Swings

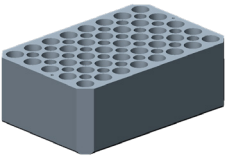
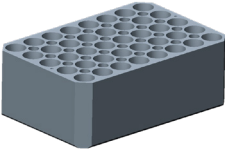
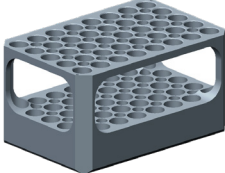
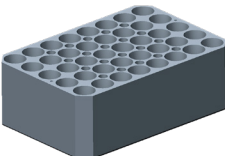
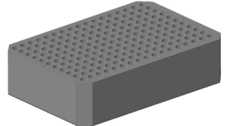
Side-bridge Swing				
	Max footprint	129.0 x 87.0 mm	Part Number	
	Mass	380 g	100022143	
	Details	Suitable for use with all systems including HCl compatible. Holds a range of sample holders.	Compatibility	
			EZ-2	✓
			S2	✓
S3			✓	

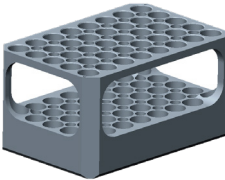
Fast-Stack Swing - Shallow Well				
	Max plate height	19 mm	Part Number	
	Max footprint	129.0 x 87.0 mm	70-1338	
	Mass	523 g	Compatibility	
	Details	Suitable for use with all systems including HCl compatible. Holds 4 shallow-well microplates.	EZ-2	✓
			S2	✓
S3			✓	

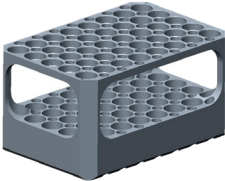
Fast-Stack Swing - Deep Well				
	Max plate height	45 mm	Part Number	
	Max footprint	129.0 x 87.0 mm	70-1612	
	Mass	400 g	Compatibility	
	Details	NOTE: Requires two heat-transfer pads (supplied). Suitable for use with all systems including HCl compatible. Holds 2 deep-well microplates. NOTE: Heat-transfer plates shall not be fitted with out microplates. NOTE: Thermal pads are consumables and be replaced when cracks are observed.	EZ-2	✓
			S2	✓
			S3	✓

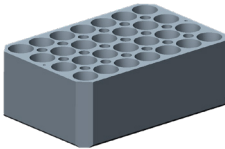
Sample Holders

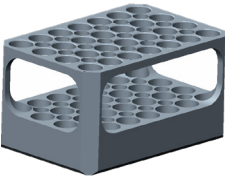
Tube holders for Side-bridge Swings

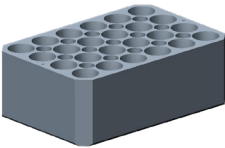
	Tube Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	10	75	10-5000	
	Limits	9.8 to 10.3	55 to 75	Compatibility	
	Details	Holds up to 40 tubes.		EZ-2	✓
				S2	✓
				S3	✓
	Tube Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	12	75	10-5025	
	Limits	11.8 to 12.3	55 to 75	Compatibility	
	Details	Holds up to 40 tubes.		EZ-2	✓
				S2	✓
				S3	✓
	Tube Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	12	100	10-5001	
	Limits	11.8 to 12.3	75 to 102	Compatibility	
	Details	Holds up to 40 tubes.		EZ-2	✓
				S2	✓
				S3	✓
	Tube Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	13	75	10-5047	
	Limits	12.8 to 13.3	55 to 75	Compatibility	
	Details	Holds up to 40 tubes.		EZ-2	✓
				S2	✓
				S3	✓
	Tube Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	6	51	10-5086	
	Limits	5.5 to 6.3	35 to 75	Compatibility	
	Details	Holds up to 187 tubes.		EZ-2	✓
				S2	✓
				S3	✓

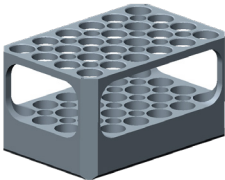
	Tube Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	13	100	10-5002	
	Limits	12.8 to 13.3	75 to 102	Compatibility	
	Details	Holds up to 40 tubes.		EZ-2	✓
				S2	✓
				S3	✓

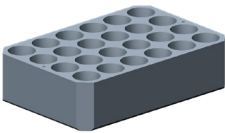
	Tube Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	13.5	100	10-5003	
	Limits	13.3 to 13.8	75 to 102	Compatibility	
	Details	Holds up to 40 tubes.		EZ-2	✓
				S2	✓
				S3	✓

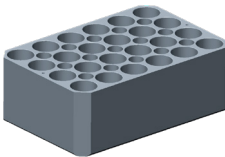
	Tube Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	15	75	10-5046	
	Limits	14.8 to 15.3	55 to 75	Compatibility	
	Details	Holds up to 28 tubes.		EZ-2	✓
				S2	✓
				S3	✓

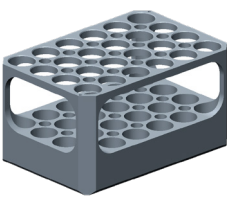
	Tube Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	15	100	10-5008	
	Limits	14.8 to 15.3	75 to 102	Compatibility	
	Details	Holds up to 28 tubes.		EZ-2	✓
				S2	✓
				S3	✓

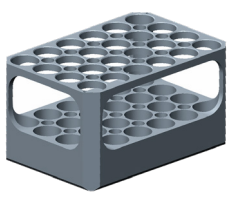
	Tube Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	16	75	10-5048	
	Limits	15.8 to 16.3	55 to 75	Compatibility	
	Details	Holds up to 24 tubes.		EZ-2	✓
				S2	✓
				S3	✓

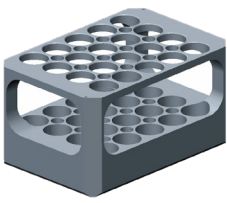
	Tube Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	16	100	10-5004	
	Limits	15.8 to 16.3	75 to 102	Compatibility	
	Details	Holds up to 24 tubes.		EZ-2	✓
				S2	✓
				S3	✓

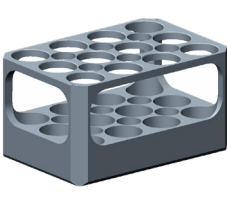
	Tube Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	16.5	50	10-5069	
	Limits	16.3 to 16.8	40 to 60	Compatibility	
	Details	Holds up to 24 tubes.		EZ-2	✓
				S2	✓
				S3	✓

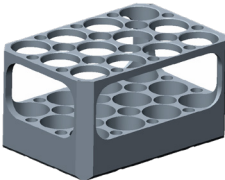
	Tube Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	17	83	10-5042	
	Limits	16.3 to 17.3	75 to 90	Compatibility	
	Details	For tapered microwave tubes. Holds up to 24 tubes.		EZ-2	✓
				S2	✓
				S3	✓

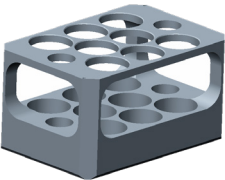
	Tube Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	17	100	10-5005	
	Limits	16.8 to 17.3	75 to 102	Compatibility	
	Details	Holds up to 24 tubes.		EZ-2	✓
				S2	✓
				S3	✓

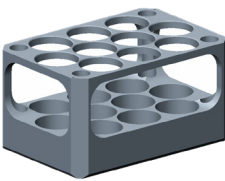
	Tube Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	18	100	10-5006	
	Limits	17.8 to 18.3	75 to 102	Compatibility	
	Details	Holds up to 24 tubes.		EZ-2	✓
				S2	✓
				S3	✓

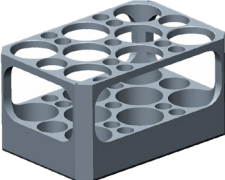
	Tube Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	18	105	10-5073	
	Limits	17.8 to 18.3	75 to 110	Compatibility	
	Details	Holds up to 20 tubes.		EZ-2	✓
				S2	✓
				S3	✓

	Tube Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	24	100	10-5044	
	Limits	23.8 to 24.3	75 to 102	Compatibility	
	Details	Holds up to 12 tubes.		EZ-2	✓
				S2	✓
				S3	✓

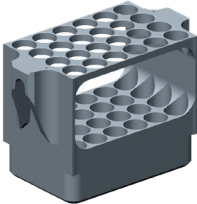
	Tube Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	25	100	10-5009	
	Limits	24.8 to 25.3	75 to 102	Compatibility	
	Details	Holds up to 12 tubes.		EZ-2	✓
				S2	✓
				S3	✓

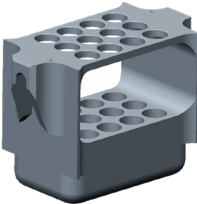
	Tube Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	27.5	100	10-5012	
	Limits	27.3 to 27.8	75 to 110	Compatibility	
	Details	Holds up to 8 tubes.		EZ-2	✓
				S2	✓
				S3	✓

	Tube Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	28	100	10-5071	
	Limits	27.8 to 28.3	75 to 110	Compatibility	
	Details	Holds up to 10 tubes.		EZ-2	✓
				S2	✓
				S3	✓

	Tube Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	30	100	10-5018	
	Limits	29.8 to 30.3	75 to 102	Compatibility	
	Details	Holds up to 8 tubes.		EZ-2	✓
				S2	✓
				S3	✓

One-Piece Tube Holders

	Tube Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	16	125	10-5070	
	Limits	15.8 to 16.3	110 to 132	Compatibility	
	Details	Holds up to 24 tubes.		EZ-2	✓
				S2	✓
				S3	✓

	Tube Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	16	150	10-5020	
	Limits	15.8 to 16.3	110 to 152	Compatibility	
	Details	Holds up to 12 tubes.		EZ-2	✓
				S2	✓
				S3	✓

	Tube Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	16.5	125	10-5050	
	Limits	16.3 to 16.8	110 to 132	Compatibility	
	Details	Holds up to 20 tubes.		EZ-2	✓
				S2	✓
				S3	✓

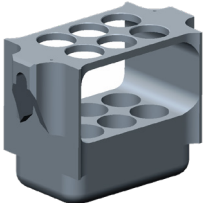
	Tube Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	17	150	10-5022	
	Limits	16.8 to 17.3	110 to 152	Compatibility	
	Details	Holds up to 24 tubes.		EZ-2	✓
				S2	✓
				S3	✓

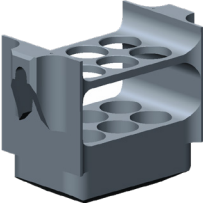
	Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	17.5	140	10-5055	
	Limits	17.0 to 17.8	105 to 142	Compatibility	
	Details	Holds up to 12 tubes.		EZ-2	✓
				S2	✓
				S3	✓

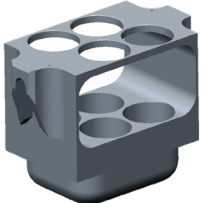
	Tube Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	18	150	10-5015	
	Limits	17.8 to 18.3	110 to 152	Compatibility	
	Details	Holds up to 12 tubes.		EZ-2	✓
				S2	✓
				S3	✓

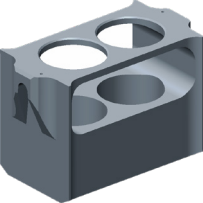
	Tube Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	20	150	10-5023	
	Limits	19.8 to 20.3	110 to 152	Compatibility	
	Details	Holds up to 10 tubes.		EZ-2	✓
				S2	✓
				S3	✓

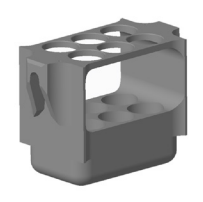
	Tube Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	24	150	10-5026	
	Limits	23.8 to 24.3	110 to 152	Compatibility	
	Details	Compatible with Radleys Carousel and STEM reaction stations. Holds up to 6 tubes.		EZ-2	✓
				S2	✓
				S3	✓

	Tube Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	25	150	10-5027	
	Limits	24.8 to 25.3	110 to 152	Compatibility	
	Details	Holds up to 6 tubes.		EZ-2	✓
				S2	✓
				S3	✓

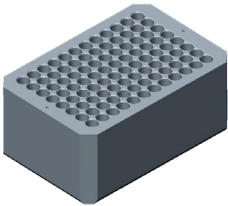
	Tube Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	26	150	10-5066	
	Limits	25.8 to 26.3	102 to 155	Compatibility	
	Details	Holds up to 6 tubes.		EZ-2	✓
				S2	✓
				S3	✓

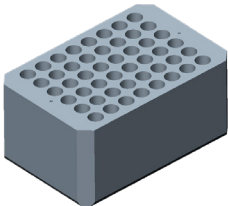
	Tube Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	34	150	10-5056	
	Limits	33.8 to 34.3	110 to 152	Compatibility	
	Details	NOTE: Requires Cushion (supplied). Holds up to 4 tubes. Replacement cushion 10-1546 (each), 10-1551 (pack of 100).		EZ-2	✓
				S2	✓
				S3	✓

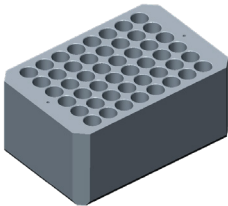
	Tube Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	45	150	10-5064	
	Limits	44.8 to 45.8	110 to 152	Compatibility	
	Details	Holds up to 2 tubes.		EZ-2	✓
				S2	✓
				S3	✓

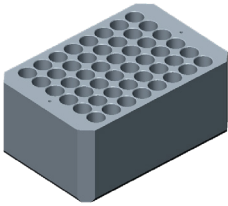
	Tube Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	27.5	140	10-5084	
	Limits	27.3 to 27.8	110 to 145	Compatibility	
	Details	Holds up to 6 tubes.		EZ-2	✓
				S2	✓
				S3	✓

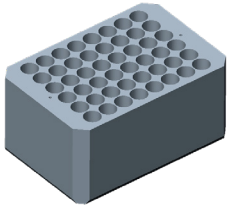
Vial-Holders for Side-Bridge Swings

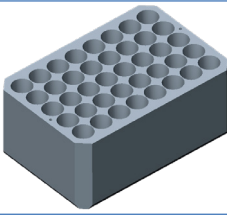
	Vial Dimensions (mm)	Diameter	Length	Part Number
	Nominal	8	35	10-5045
	Limits	7.5 to 8.3	30 to 40	Compatibility
	Details	Holds up to 96 vials.		EZ-2 ✓
				S2 ✓
				S3 ✓

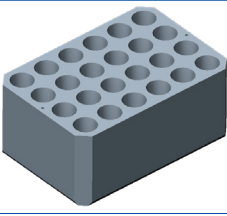
	Vial Dimensions (mm)	Diameter	Length	Part Number
	Nominal	10	45	10-5029
	Limits	9.5 to 10.3	35 to 45	Compatibility
	Details	Holds up to 48 vials.		EZ-2 ✓
				S2 ✓
				S3 ✓

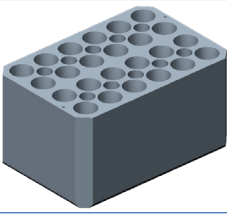
	Vial Dimensions (mm)	Diameter	Length	Part Number
	Nominal	11	35	10-5032
	Limits	10.7 to 11.5	30 to 40	Compatibility
	Details	Holds up to 48 vials.		EZ-2 ✓
				S2 ✓
				S3 ✓

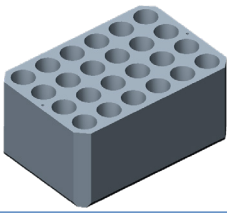
	Vial Dimensions (mm)	Diameter	Length	Part Number
	Nominal	11.5	55	10-5074
	Limits	11.0 to 11.8	40 to 60	Compatibility
	Details	Holds up to 48 vials.		EZ-2 ✓
				S2 ✓
				S3 ✓

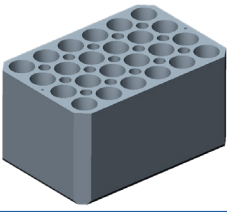
	Vial Dimensions (mm)	Diameter	Length	Part Number
	Nominal	12	35	10-5024
	Limits	11.3 to 12.1	30 to 40	Compatibility
	Details	Including GC auto-sampler vials. Holds up to 48 vials.		EZ-2 ✓
				S2 ✓
				S3 ✓

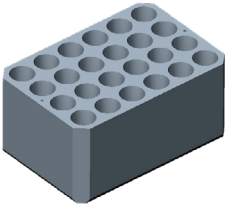
	Vial Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	13	50	10-5034	
	Limits	12.5 to 13.3	30 to 50	Compatibility	
	Details	Holds up to 40 vials.		EZ-2	✓
				S2	✓
				S3	✓

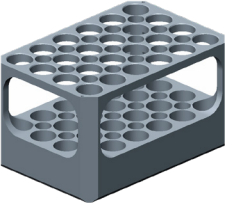
	Vial Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	14.5	45	10-5035	
	Limits	14.0 to 14.8	35 to 50	Compatibility	
	Details	Holds up to 24 vials.		EZ-2	✓
				S2	✓
				S3	✓

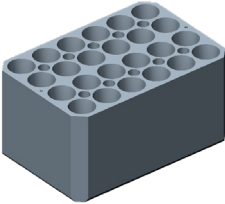
	Vial Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	14.5	65	10-5016	
	Limits	14.0 to 14.8	55 to 75	Compatibility	
	Details	Holds up to 24 vials.		EZ-2	✓
				S2	✓
				S3	✓

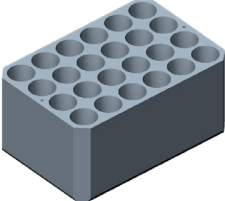
	Vial Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	15	45	10-5037	
	Limits	14.3 to 15.1	35 to 50	Compatibility	
	Details	Holds up to 24 vials.		EZ-2	✓
				S2	✓
				S3	✓

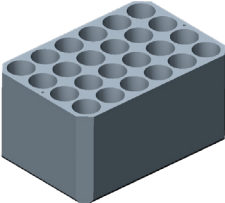
	Vial Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	15	65	10-5021	
	Limits	14.3 to 15.1	55 to 75	Compatibility	
	Details	Holds up to 24 vials.		EZ-2	✓
				S2	✓
				S3	✓

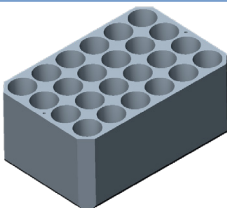
	Vial Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	15.5	45	10-5038	
	Limits	15.0 to 15.8	35 to 50	Compatibility	
	Details	Holds up to 24 vials.		EZ-2	✓
				S2	✓
				S3	✓

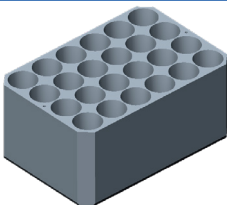
	Vial Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	16	100	10-5067	
	Limits	15.5 to 16.3	75 to 102	Compatibility	
	Details	Holds up to 24 vials.		EZ-2	✓
				S2	✓
				S3	✓

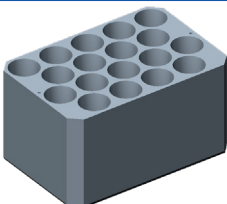
	Vial Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	16.5	65	10-5028	
	Limits	16.2 to 17.0	55 to 75	Compatibility	
	Details	Holds up to 24 vials.		EZ-2	✓
				S2	✓
				S3	✓

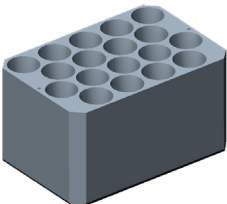
	Vial Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	17	40	10-5007	
	Limits	16.3 to 17.1	38 to 50	Compatibility	
	Details	Holds up to 24 vials.		EZ-2	✓
				S2	✓
				S3	✓

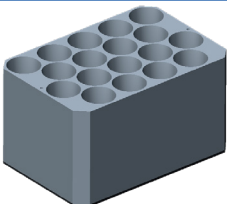
	Vial Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	17	65	10-5030	
	Limits	16.5 to 17.3	55 to 75	Compatibility	
	Details	Including 2 dram vial. Holds up to 24 vials.		EZ-2	✓
				S2	✓
				S3	✓

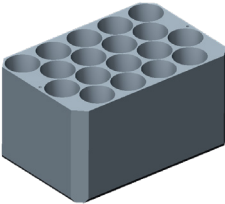
	Vial Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	17.5	45	10-5031	
	Limits	17.0 to 17.8	40 to 50	Compatibility	
	Details	Holds up to 24 vials.		EZ-2	✓
				S2	✓
				S3	✓

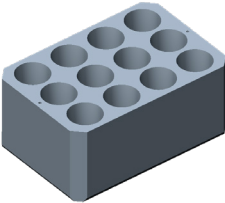
	Vial Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	18	45	10-5033	
	Limits	17.7 to 18.5	40 to 50	Compatibility	
	Details	Holds up to 24 vials.		EZ-2	✓
				S2	✓
				S3	✓

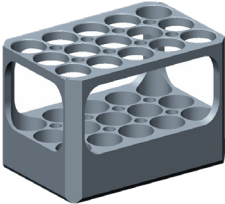
	Vial Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	20	65	10-5036	
	Limits	19.5 to 20.3	55 to 75	Compatibility	
	Details	Holds up to 18 vials.		EZ-2	✓
				S2	✓
				S3	✓

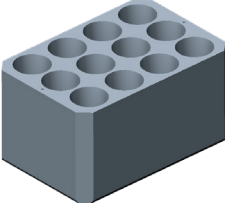
	Vial Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	20.5	65	10-5040	
	Limits	20.0 to 20.8	55 to 75	Compatibility	
	Details	Holds up to 18 vials.		EZ-2	✓
				S2	✓
				S3	✓

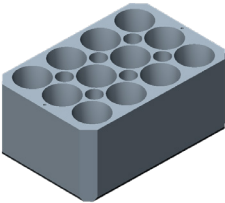
	Vial Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	21	65	10-5039	
	Limits	20.5 to 21.3	55 to 75	Compatibility	
	Details	Holds up to 18 vials.		EZ-2	✓
				S2	✓
				S3	✓

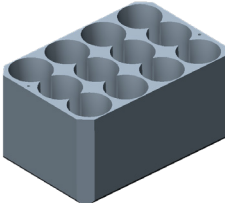
	Vial Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	21.5	65	10-5041	
	Limits	21.0 to 21.8	55 to 75	Compatibility	
	Details	Holds up to 18 vials.		EZ-2	✓
				S2	✓
				S3	✓

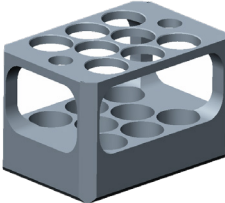
	Vial Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	22.5	50	10-5062	
	Limits	22.0 to 22.8	38 to 50	Compatibility	
	Details	Holds up to 12 vials.		EZ-2	✓
				S2	✓
				S3	✓

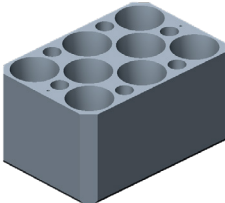
	Vial Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	23	85	10-5079	
	Limits	22.5 to 23.3	75 to 102	Compatibility	
	Details	Holds up to 15 vials.		EZ-2	✓
				S2	✓
				S3	✓

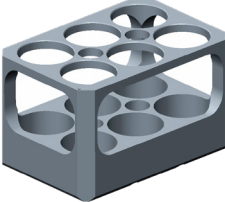
	Vial Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	23.5	65	10-5019	
	Limits	23.0 to 23.8	55 to 75	Compatibility	
	Details	Holds up to 12 vials.		EZ-2	✓
				S2	✓
				S3	✓

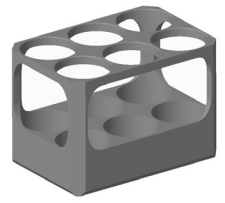
	Vial Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	24	45	10-5075	
	Limits	23.7 to 24.5	45 to 75	Compatibility	
	Details	Holds up to 12 vials.		EZ-2	✓
				S2	✓
				S3	✓

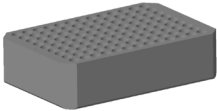
	Vial Dimensions (mm)	Diameter	Length	Part Number
	Nominal	27.5	65	10-5013
	Limits	27.0 to 27.8	55 to 75	Compatibility
	Details	Accepts 20 ml scintillation vials. Holds up to 12 vials.		EZ-2 ✓
				S2 ✓
				S3 ✓

	Vial Dimensions (mm)	Diameter	Length	Part Number
	Nominal	27.5	100	10-5010
	Limits	27.0 to 27.8	75 to 110	Compatibility
	Details	NOTE: Requires Cushion (supplied). Accepts 40 ml scintillation vials. Holds up to 8 vials. Replacement cushion 10-1546 (each), 10-1551 (pack of 100).		EZ-2 ✓
				S2 ✓
				S3 ✓

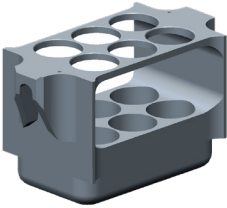
	Vial Dimensions (mm)	Diameter	Length	Part Number
	Nominal	30	65	10-5078
	Limits	29.5 to 30.3	55 to 75	Compatibility
	Details	Holds up to 8 vials.		EZ-2 ✓
				S2 ✓
				S3 ✓

	Vial Dimensions (mm)	Diameter	Length	Part Number
	Nominal	35	100	10-5068
	Limits	34.5 to 35.8	75 to 102	Compatibility
	Details	Holds up to 6 vials.		EZ-2 ✓
				S2 ✓
				S3 ✓

	Vial Dimensions (mm)	Diameter	Length	Part Number
	Nominal	40	75	10-5085
	Limits	39.5 to 40.4	72 to 80	Compatibility
	Details	Holds up to 6 vials.		EZ-2 ✓
				S2 ✓
				S3 ✓

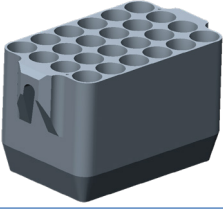
	Vial Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	7	32	10-5087	
	Limits	6.5 to 7.5	22 to 42	Compatibility	
	Details	Holds up to 117 vials.		EZ-2	✓
				S2	✓
				S3	✓

One-Piece Vial Holders

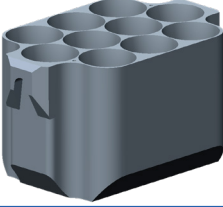
	Vial Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	27.5	140	10-5076	
	Limits	27.0 to 27.8	105 to 142	Compatibility	
	Details	NOTE: Requires thermal cushion (10-1546). For Dionex 60 ml ASE® tubes. Holds up to 6 vials. Replacement cushion 10-1546 (each), 10-1551 (pack of 100).		EZ-2	✓
				S2	✓
				S3	✓

ASE® is a registered trademark of Dionex Corporation.

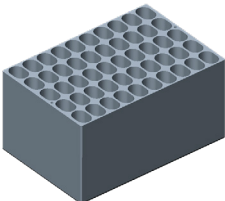
Centrifuge Tube Holders

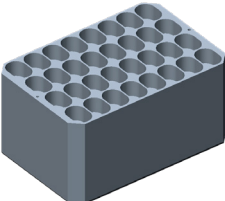
	Tube Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	16.5	115	70-1192	
	Limits	16.3 to 17.3	100 to 120	Compatibility	
	Details	NOTE: Requires tip cushion (supplied), Holds up to 24 tubes. Replaces 10-5053.		EZ-2	✓
				S2	✓
				S3	✓

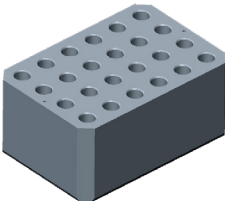
For 30° ended 15 ml centrifuge tubes.
 Replacement cushion 10-1585 (each), 10-1602 (pack of 24).

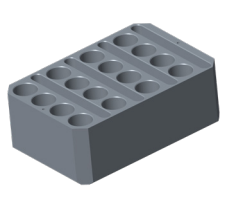
	Tube Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	29.5	115	10-5054	
	Limits	28.5 to 29.8	100 to 115	Compatibility	
	Details	For 70° ended 50 ml centrifuge tubes. Holds up to 10 tubes.		EZ-2	✓
				S2	✓
				S3	✓

Microcentrifuge Tube Holders

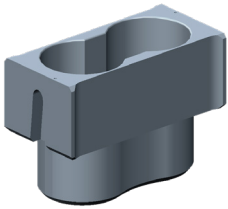
	Tube Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	8	30	10-5049	
	Limits	NA	NA	Compatibility	
	Details	For 0.5 ml flip-cap, without skirt. Holds up to 50 tubes.		EZ-2	✓
				S2	✓
				S3	✓

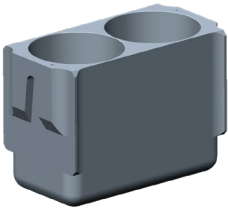
	Tube Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	10	40	10-5043	
	Limits	NA	NA	Compatibility	
	Details	For 1.5 ml flip-cap, without skirt. Holds up to 32 tubes.		EZ-2	✓
				S2	✓
				S3	✓

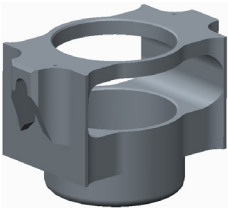
	Tube Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	8	30	10-5011	
	Limits	NA	NA	Compatibility	
	Details	For 2.0 ml flip-cap, without skirt. Holds up to 24 tubes.		EZ-2	✓
				S2	✓
				S3	✓

	Tube Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	17	40	10-5089	
	Limits	NA	NA	Compatibility	
	Details	For 5.0 ml flip-cap, without skirt. Holds up to 16 tubes.		EZ-2	✓
				S2	✓
				S3	✓

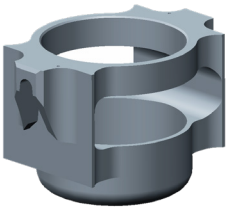
Bottle Holders

	Bottle Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	55.5	100	70-1025	
	Limits	54.0 to 56.0	90 to 110	Compatibility	
	Details	Including Schott 100 ml. Holds 2 bottles.		EZ-2	✓
				S2	✓
				S3	✓

	Bottle Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	60.5	125	10-5061	
	Limits	59.5 to 62	105 to 127	Compatibility	
	Details	Holds 2 bottles.		EZ-2	✓
				S2	✓
				S3	✓

	Bottle Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	70	140	70-1795	
	Limits	69 to 70.8	110 to 140	Compatibility	
	Details	NOTE: Requires cushion (supplied). Replacement cushion: 04-8963.		EZ-2	✓
				S2	✓
				S3	✓

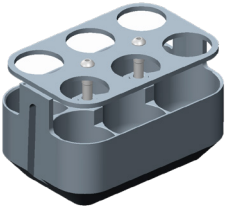
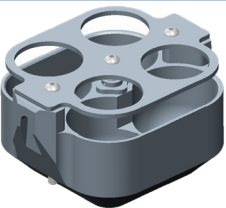
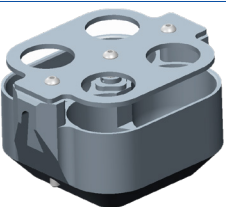
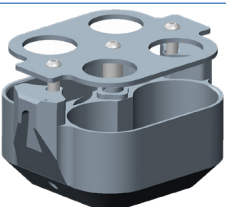
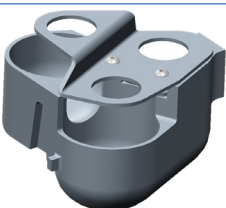
Including Schott 250ml & 400ml beaker.

	Bottle Dimensions (mm)	Diameter	Length	Part Number	
	Nominal	82	152	10-5063	
	Limits	81 to 82.8	110 to 152	Compatibility	
	Details	NOTE: Not for use with EZ-2 Mark 1 models. Requires cushion (supplied).		EZ-2	✓
				S2	✓
				S3	✓

For 500 ml bottle. Holds one bottle.
Replacement cushion 10-1543 (each), 10-1553 (pack of 100).

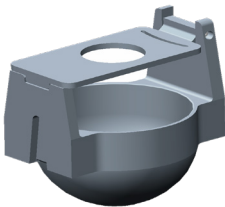
Flask Holders

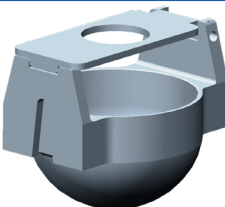
Flask socket sizes are European ISO-standard joint sizes. The first number represents the outer diameter at the widest part of the ground glass joint section in millimetres. The second number represents the length of the joint section in millimetres.

	Flask Volume	25 ml	Part Number	
	Socket Size	19/26 to 24/29	70-0538	
	Details	NOTE: Requires cushion (supplied). Holds 6 flasks. Replacement cushion 10-1546 (each), 10-1551 (pack of 100).	Compatibility	
			EZ-2	✓
			S2	✓
			S3	✓
	Flask Volume	50 ml	Part Number	
	Socket Size	29/32	70-0539	
	Details	NOTE: Requires cushion (supplied). Holds 4 flasks. Replacement cushion 10-1540 (each), 10-1556 (pack of 100).	Compatibility	
			EZ-2	✓
			S2	✓
			S3	✓
	Flask Volume	50 ml	Part Number	
	Socket Size	19/36 to 24/29	70-0656	
	Details	NOTE: Requires cushion (supplied). Holds 4 flasks. Replacement cushion 10-1540 (each), 10-1556 (pack of 100).	Compatibility	
			EZ-2	✓
			S2	✓
			S3	✓
	Flask Volume	50 ml	Part Number	
	Socket Size	24/29 to 29/32	70-1026	
	Details	NOTE: Requires cushion (supplied). Holds 4 flasks. Replacement cushion 10-1540 (each), 10-1556 (pack of 100).	Compatibility	
			EZ-2	✓
			S2	✓
			S3	✓
	Flask Volume	100 ml	Part Number	
	Socket Size	19/26 to 24/29	70-0657	
	Details	NOTE: Requires cushion (supplied). Compatible with Radleys 6-piece Carousel. Holds 3 flasks. Replacement cushion 10-1541 (each), 10-1555 (pack of 48).	Compatibility	
			EZ-2	✓
			S2	✗
			S3	✗

	Flask Volume	100 ml	Part Number
	Socket Size	29/36 to 19/32	70-0579
	Details	NOTE: Requires cushion (supplied). Holds 3 flasks. Replacement cushion 10-1541 (each), 10-1555 (pack of 48).	Compatibility
			EZ-2 ✓
			S2 ✗
			S3 ✗

	Flask Volume	250 ml	Part Number
	Socket Size	19/22 to 29/32	70-0580
	Details	NOTE: Requires cushion (supplied). Compatible with Radleys 6-piece Carousel. Holds 1 flask. Replacement cushion 10-1543 (each), 10-1553 (pack of 12).	Compatibility
			EZ-2 ✓
			S2 ✓
			S3 ✓

	Flask Volume	500 ml	Part Number
	Socket Size	19/32 to 29/32	70-0582
	Details	NOTE: Requires cushion (supplied). Holds 1 flask. Replacement cushion 10-1544 (each), 10-1552 (pack of 12).	Compatibility
			EZ-2 ✓
			S2 ✗
			S3 ✗

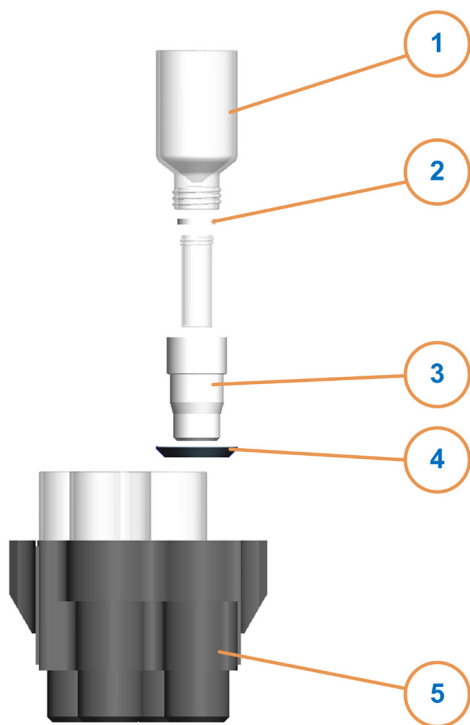
	Flask Volume	500 ml	Part Number
	Socket Size	19/32 to 29/32	10-1616
	Details	NOTE: Requires cushion (supplied). Max. flask height: 156mm. Holds 1 flask. Replacement cushion 10-1544 (each), 10-1552 (pack of 12).	Compatibility
			EZ-2 ✗
			S2 ✓
			S3 ✓

SampleGenie™

The revolutionary SampleGenie technology enables large volume samples to be dried or concentrated directly into a variety of small vials. Where HPLC purification fractions are used, fractions are combined into one tube and dried or fast lyophilised into the final vial.



60 ml SampleGenie



Item	Description	Qty in Set	Part Number
1	Flask	10	70-1153
2	Seal for GC vials - white Viton	▪	70-1359
	Seal for GC vials - perfluoroelastomer	▪	70-1360
	Seal for all vials except GC - white Viton	▪	70-1361
	Seal for all vials except GC - perfluoroelastomer	▪	70-1362
3	Adapter	▪	▪ ▪
4	Cushion for 60 ml SampleGenie - Viton	10	70-1448
5	Sample Holder:		
	SampleGenie set with white Viton seals	1	70-1340
	SampleGenie set with perfluoroelastomer seals	1	70-1341

▪ Specify quantity when ordering

▪ ▪ Refer to following table for 60 ml adapter part numbers

Please specify adapter type when ordering complete assemblies.

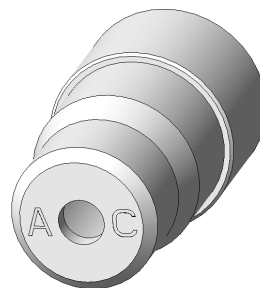
60 ml Drying Adapters

Code	Max. Vial Diameter (mm)	Vial Length (mm)	Part Number
AE	11.8	30.5 to 33.5	04-6214
AG	11.8	45.5 to 48.5	04-6615
AD	13.5	38.5 to 41.5	04-6174
AL	15.3	30.5 to 33.5	04-7593
AB	15.3	43.5 to 46.5	04-6155
AF	15.3	48.5 to 51.5	04-4686
AK	15.6	38.5 to 41.5	04-7302
AM	16.6	48.5 to 51.5	15-0448
AJ	16.8	43.5 to 46.5	04-6756
AA	17.3	48.5 to 51.5	04-6156
AH	19.8	38.5 to 41.5	04-6619
AC	20.3	48.5 to 51.1	04-6166



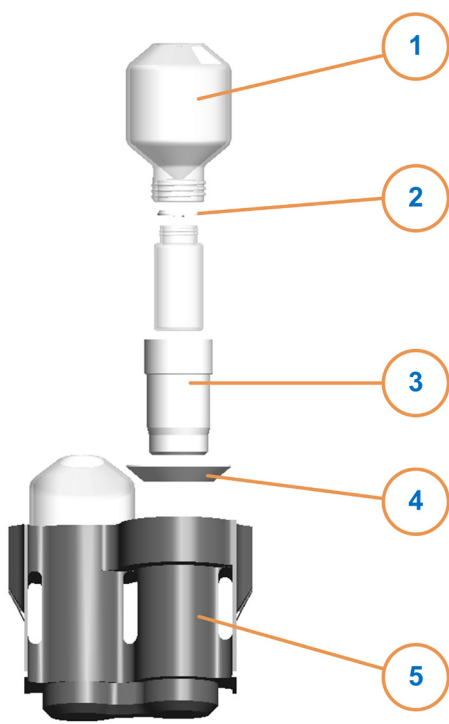
60 ml Lyophilisation/Concentration Adapters

Code	Max. Vial Diameter (mm)	Vial Length (mm)	Part Number
AE	11.8	30.5 to 33.5	04-6215
AG	11.8	45.5 to 48.5	04-6614
AD	13.5	38.5 to 41.5	04-6175
AL	15.3	30.5 to 33.5	04-7592
AB	15.3	43.5 to 46.5	04-6130
AF	15.3	48.5 to 51.5	04-4688
AK	15.6	38.5 to 41.5	04-7301
AM	16.6	48.5 to 51.5	15-0450
AJ	16.8	43.5 to 46.5	04-6757
AA	17.3	48.5 to 51.5	04-6110
AH	19.8	38.5 to 41.5	04-6618
AC	20.3	48.5 to 51.1	04-6167



NOTE: Code is shown on the base of the adaptor.

125 ml SampleGenie



Item	Description	Qty in Set	Part Number
1	Flask	10	70-1157
2	Seal for GC vials - white Viton	▪	70-1359
	Seal for GC vials - perfluoroelastomer	▪	70-1360
	Seal for all vials except GC - white Viton	▪	70-1363
	Seal for all vials except GC - perfluoroelastomer	▪	70-1364
3	Adapter	▪	▪ ▪
4	Cushion for 125 ml SampleGenie - PTFE	8	70-1373
5	Sample Holder:		
	SampleGenie set with white Viton seals	1	70-1346
	SampleGenie set with perfluoroelastomer seals	1	70-1347

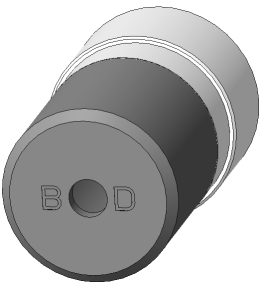
▪ Specify quantity when ordering

▪ ▪ Refer to following table for 125 ml adapter part numbers

Please specify adapter type when ordering complete assemblies.

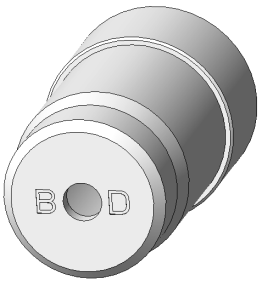
125 ml Drying Adapters

Code	Max. Vial Diameter (mm)	Vial Length (mm)	Part Number
BE	11.8	30.5 to 33.5	04-6216
BC	15.3	43.5 to 46.5	04-6162
BG	15.3	48.5 to 51.5	04-6378
BB	16.8	33.5 to 36.5	04-6161
BM	16.8	43.5 to 46.5	04-6759
BH	17.3	48.5 to 51.5	04-6390
BF	17.3	58.5 to 61.5	04-6292
BL	18.0	43.5 to 46.5	04-6737
BO	18.3	38.5 to 41.5	04-7645
BK	21.8	55.5 to 58.5	04-6606
BJ	23.8	58.5 to 61.5	04-6454
BN	24.3	48.5 to 51.5	04-6979
BA	27.3	58.5 to 61.5	04-6048
BD	28.3	58.5 to 61.5	04-6190



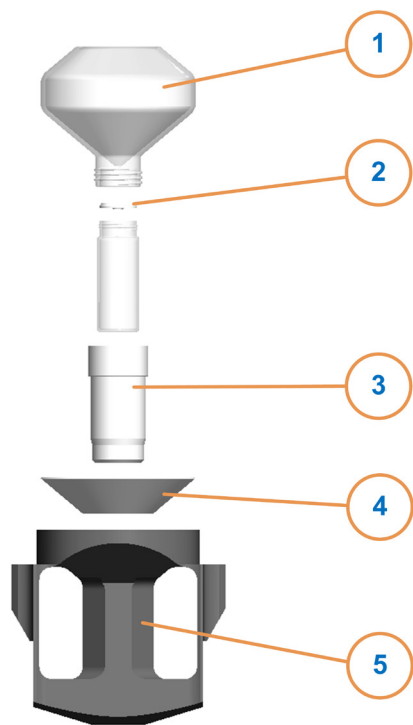
125 ml Lyophilisation/Concentration Adapters

Code	Max. Vial Diameter (mm)	Vial Length (mm)	Part Number
BE	11.8	30.5 to 33.5	04-6217
BC	15.3	43.5 to 46.5	04-6160
BG	15.3	48.5 to 51.5	04-6377
BB	16.8	33.5 to 36.5	04-6159
BM	16.8	43.5 to 46.5	04-6760
BH	17.3	48.5 to 51.5	04-6389
BF	17.3	58.5 to 61.5	04-6293
BL	18.0	43.5 to 46.5	04-6739
BO	18.3	38.5 to 41.5	04-7644
BK	21.8	55.5 to 58.5	04-6605
BJ	23.8	58.5 to 61.5	04-6453
BN	24.3	48.5 to 51.5	04-6980
BA	27.3	58.5 to 61.5	04-6049
BD	28.3	58.5 to 61.5	04-6192



NOTE: Code is shown on the base of the adaptor.

250 ml SampleGenie



Item	Description	Qty in Set	Part Number
1	Flask	4	70-1158
2	Seal for GC vials - white Viton	▪	70-1359
	Seal for GC vials - perfluoroelastomer	▪	70-1360
	Seal for all vials except GC - white Viton	▪	70-1363
	Seal for all vials except GC - perfluoroelastomer	▪	70-1364
3	Adapter	▪	▪ ▪
4	Cushion for 125 ml SampleGenie - PTFE	4	70-1358
5	Sample Holder:		
	SampleGenie set with white Viton seals	1	70-1352
	SampleGenie set with perfluoroelastomer seals	1	70-1353

▪ Specify quantity when ordering

▪ ▪ Refer to following table for 250 ml adapter part numbers

Please specify adapter type when ordering complete assemblies.

250 ml Drying Adapters

Code	Max. Vial Diameter (mm)	Vial Length (mm)	Part Number
CF	11.8	30.5 to 33.5	04-6218
CE	15.3	43.5 to 46.5	04-6186
CD	17.3	58.5 to 61.5	04-6184
CH	21.3	68.5 to 71.5	04-6363
CC	27.3	58.5 to 61.5	04-6182
CA	27.3	68.5 to 71.5	04-6163
CG	28.3	58.5 to 61.5	04-6238
CB	28.3	68.5 to 71.5	04-6168



250 ml Lyophilisation/Concentration Adapters

Code	Max. Vial Diameter (mm)	Vial Length (mm)	Part Number
CF	11.8	30.5 to 33.5	04-6219
CE	15.3	43.5 to 46.5	04-6172
CD	17.3	58.5 to 61.5	04-6171
CH	21.3	68.5 to 71.5	04-6245
CC	27.3	58.5 to 61.5	04-6170
CA	27.3	68.5 to 71.5	04-6144
CG	28.3	58.5 to 61.5	04-6239
CB	28.3	68.5 to 71.5	04-6169



NOTE: Code is shown on the base of the adaptor.

Miscellaneous

Infinity Trolley

The Infinity Trolley is a modular system designed to accommodate a range of SP Genevac evaporation equipment. The trolley provides safe and durable support for the evaporator, whilst occupying the minimum amount of laboratory space. The lower “truck” can be extended to provide access to items on the lower shelf.



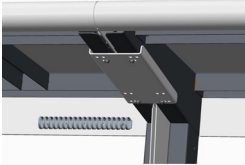
Approximate Infinity Trolley Dimensions:

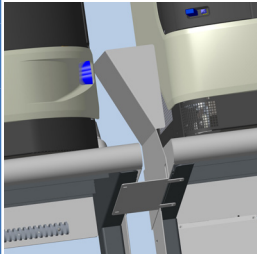
740 mm x 733 mm x 835 mm (W x D x H)

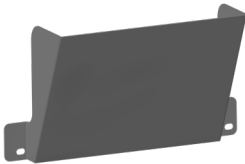
When truck is fully extended, add 500 mm to depth.
For optional side-shelf, add 260 mm to width.

	Infinity Trolley		Part Number
	Description	Infinity Trolley	70-1850
	Details	Compatible with miVac, EZ-2 Series, Rocket Synergy and Series 3 evaporators.	

	Infinity Trolley Accessory		Part Number
	Description	Shelf Kit	70-1635
	Details	Bolts to the back of the Infinity Trolley. May be fitted to left- and right-hand sides.	

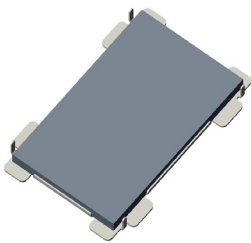
	Infinity Trolley Accessory		Part Number
	Description	Joining Kit	70-1636
	Details	Plate and fixing screws to securely join an Infinity Trolley to another adjacent Trolley. Any number of Infinity Trollies may be joined together, using one joining kit for each Trolley being joined. NOTE: Not suitable for use with Series 3 systems.	

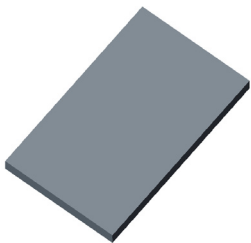
	Infinity Trolley Accessory		Part Number
	Description	Series 3 Deflector Plate & Joining Kit	70-1900
	Details	Designed specifically for the Series 3, this Infinity Trolley joining kit includes a deflector plate to direct exhaust airflow away from the evaporator. NOTE: Suitable for use with Series 3 systems only.	

	Series 3 Accessory		Part Number
	Description	Series 3 Deflector Plate	70-2017
	Details	For use with Series 3 evaporators. Directs exhaust airflow away from adjacent evaporators when siting multiple Series 3 systems together on a bench (without Infinity Trollies).	

Heat-Transfer Plates for Microplates

Heat-Transfer Plates provide improved evaporation results when using microtitre plates. Heat-Transfer Plates are suitable for use with microtitre plates only.

	Heat-Transfer Plates		Part Number
	Description	HTP - Complete Assembly	70-1786
	Details	Heat-Transfer Plate with pliable thermal pad. Compatible with all types of deep-well plate that fit within the footprint of the fast-stack or side-bridge swing.	

	Heat-Transfer Plates		Part Number
	Description	HTP - Thermal Pad	04-8824
	Details	Replacement thermal pad (only) for Heat-Transfer Plate Assembly.	

NOTE: Heat-Transfer Plate assembly 70-1786 is supplied with deep-well fast-stack swings (70-1612), and may also be used with standard side-bridge swings (100022143).

NOTE: Thermal Pads are a consumable item. They may be turned over, allowing both sides of the pad to be used before replacement.

Highly durable, solid aluminium Heat-Transfer Plates are available to suit a range of microtitre plates. Please contact your local Genevac representative for more details.



NOTE: Solid Heat-Transfer Plates should only be used with the type of microtitre plate for which they were specified.

Gilson and Waters Fraction Collection Racks

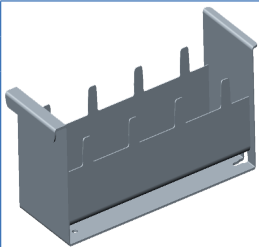
Genevac offer a range of innovative options that help interface fraction collection systems directly to a Genevac evaporator. Gilson 215 and Waters 2767 systems are currently supported.

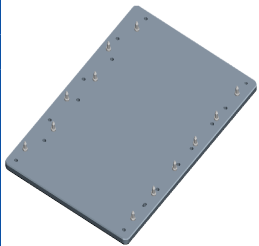
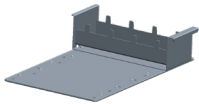
To interface with the fraction collection system, SP provide a simple sample holder carrier which fits directly onto the fraction collection bed. SP Genevac sample holders fit into the carrier to make up a complete fraction collection rack. Once fractions are collected, the sample holders can be placed directly into the SP Genevac evaporator with no need to label, mark or manipulate individual tubes. Sample holders can be marked to indicate their location within the fraction collection rack.

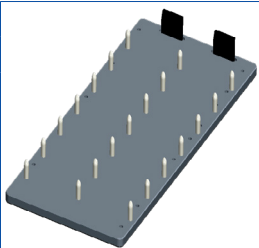
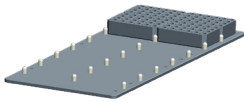
A range of sample holders is available, including standard sample holders and specially-designed holders which ensure uniform spacing between tubes and help reduce fraction collector re-programming.

NOTE: Waters 2767 users will require the adapter plate when using 16 mm diameter tubes.

Further information, including a technical article discussing Fraction Collection Racks in greater detail, is available upon request from your local SP representative.

	Rack		Part Number
	Description	Rack	10-1570
	Details	For use with Waters Fraction Collection Rack. Holds 4 tube holders (e.g. 10-5020 or 10-5027).	

	Adapter Plate		Part Number
	Description	Adapter Plate for 10-1570	04-4181
	Details	Converts Waters 2767 Sample Manager Rack Bed to hold 3 racks. 	

	Adapter Plate		Part Number
	Description	Plate for side-bridge holders	04-4330
	Details	Converts Waters 2767 Sample Manager Rack Bed to hold 14 MAX side-bridge holders. 	

NOTE: The items listed above represent a small sample from the range of fraction collection racks that Genevac can provide. Please contact your local Genevac representative for more details.

Thermal Cushions

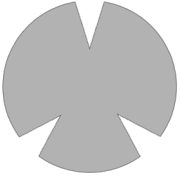
Genevac Thermal Cushions prolong glassware lifetime and dramatically reduce sample drying time. The use of Thermal Cushions is mandatory with some sample holders.

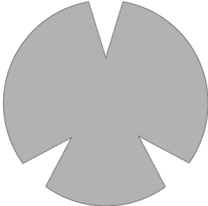
	Thermal Cushion		Part Number
	Diameter	25 mm	10-1565
	Details	Special. Pack of 100.	

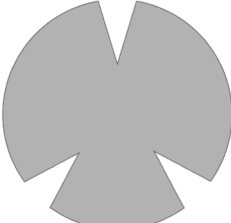
	Thermal Cushion		Part Number
	Diameter	28 mm	10-1551
	Details	For use with flasks or tubes up to 25 ml. Pack of 100.	

	Thermal Cushion		Part Number
	Diameter	33 mm	10-1556
	Details	For use with flasks up to 50 ml. Pack of 100.	

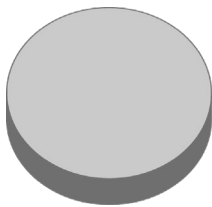
	Thermal Cushion		Part Number
	Diameter	38 mm	10-1554
	Details	For use with flasks up to 170 ml. Pack of 48.	

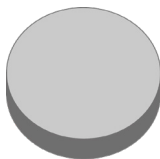
	Thermal Cushion		Part Number
	Diameter	42 mm	10-1555
	Details	For use with flasks up to 100 ml. Pack of 48.	

	Thermal Cushion		Part Number
	Diameter	68 mm	10-1553
	Details	For use with flasks up to 250 ml. Pack of 12.	

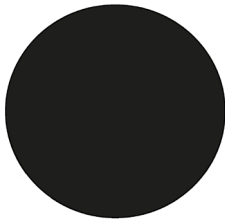
	Thermal Cushion		Part Number
	Diameter	82 mm	10-1552
	Details	For use with flasks up to 500 ml. Pack of 12.	

Tip Cushions

	Tip Cushion		Part Number
	Quantity	Set of 24	70-1371
	Details	For use with sample holder 10-5065.	

	Tip Cushion		Part Number
	Quantity	Set of 24	10-1602
	Details	For use with sample holder 70-1004.	

Viton Cushion

	Viton Thermal Cushion		Part Number
	Diameter	68 mm	04-8963
	Details	For use with sample holders 70-1795 and 10-5060. Please specify quantity required when ordering.	

Spare Parts

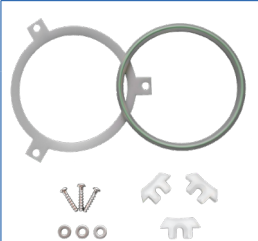
EZ-2 Consumables

The following parts are compatible with all EZ-2 series evaporators that have a detachable condenser jar.

EZ-2 Mk2 & Mk3

	Spare Part		Part Number
	Description	Infra-Red Lamp	AC9060
	Details	For use with EZ-2.2, EZ-2.3 and EZ-2 Elite. Lamp only.	

	Spare Part		Part Number
	Description	Lid Seal	04-5094/S
	Details	For use with EZ-2.2, EZ-2.3 and EZ-2 Elite.	

	Spare Part		Part Number
	Description	Jar Seal Kit	70-1494/S
	Details	For use with EZ-2.2 and EZ-2.3. Seal and thread insert kit for condenser jar.	

	Spare Part		Part Number
	Description	Condenser Jar	STJ-10000-000
	Details	For use with EZ-2.2 and EZ-2.3. Spare condenser jar.	

NOTE: For fitting instructions, please refer to the EZ-2 User Manual.

EZ-2 4.0

	Spare Part		Part Number
	Description	High Power Infra-Red Lamp	100022144
	Details	For use with EZ-2 4.0. Lamp only.	

	Spare Part		Part Number
	Description	Lid Seal	100019000
	Details	For use with EZ-2 4.0.	

	Spare Part		Part Number
	Description	Condenser Seal	100018831
	Details	For use with EZ-2 4.0.	

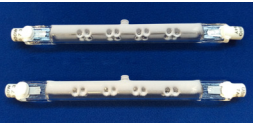
	Spare Part		Part Number
	Description	Condenser Jar	100022176
	Details	For use with EZ-2 4.0 Standard & Plus. Spare condenser jar.	

	Spare Part		Part Number
	Description	Waste Container, 3-Port	70-1656/S
	Details	For use with EZ-2 4.0 Elite.	

	Spare Part		Part Number
	Description	Waste Container, 2-Port	70-1386/S
	Details	For use with EZ-2 4.0 Elite with Inert Gas Purge Option.	

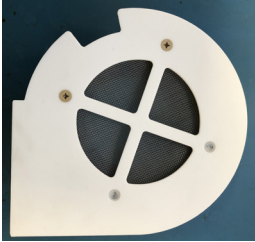
Series 3 Consumables

	Series 3 Chamber Door Seal		Part Number
	Dimensions	435 mm x 444 mm (W x H)	70-2050/S
	Details	For Series 3 evaporators.	Replacement chamber door seal for Series 3 evaporators.

	Series 3 Infra-Red Lamps		Part Number
	Quantity	Set of two.	70-1846/S
	Details	For Series 3 evaporators.	Replacement IR lamps for Series 3 evaporators.

	Series 3 SampleGuard Battery Housing Seal		Part Number
	Quantity	Set of one.	70-1848/S
	Details	For Series 3 evaporators.	Replacement SampleGuard battery housing seal for Series 3 evaporators.

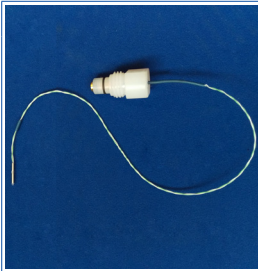
	Series 3 SampleGuard Probe Sealing O-Ring		Part Number
	Quantity	Set of four.	70-1847/S
	Details	For Series 3 evaporators.	Replacement SampleGuard probe sealing o-rings for Series 3 evaporators.

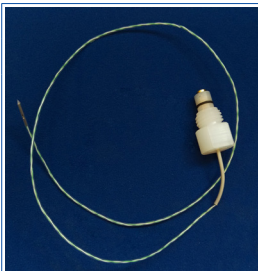
	Series 3 Condenser Filter		Part Number
	Quantity	Each.	70-1851/S
	Details	For Series 3 evaporators.	Replacement condenser filter for Series 3 evaporators.

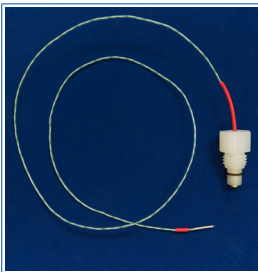
NOTE: For fitting instructions, please refer to the *Series 3 Installation & Maintenance Manual*.

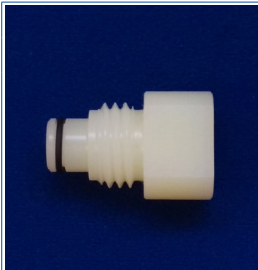
Series 3 SampleGuard Probes

All Series 3 SampleGuard probes are HCl-compatible. Probes are a consumable item.

	Series 3 SampleGuard Probes		Part Number
	Length	445 mm	70-2091/S
	Details	For Series 3 evaporators.	Colour: Blue For Rotor position A1 (Control)

	Series 3 SampleGuard Probes		Part Number
	Length	655 mm	70-2090/S
	Details	For Series 3 evaporators.	Colour: Grey For Rotor positions A1 (Auxiliary) and B4 (Control)

	Series 3 SampleGuard Probes		Part Number
	Length	765 mm	70-2092/S
	Details	For Series 3 evaporators.	Colour: Red For Rotor position B4 (Auxiliary)

	Series 3 SampleGuard Probe Blanking Plug		Part Number
	Quantity		70-1852/S
	Details	For Series 3 evaporators.	Replacement SampleGuard Probe Blanking Plug for Series 3 evaporators.

NOTE: For fitting instructions, please refer to the *Series 3 Installation & Maintenance Manual*.

This page intentionally left blank.

This page intentionally left blank.

This page intentionally left blank.

Genevac Patents

Genevac products are protected by the following patents and patent applications.

1297873	2332865	1297873 FR	2396575
1153278	2334688	1556149 FR	2363745
1073505 FR	6,605,474	1556149	2332865
699 18 734.6-08	1297874	600 07 636.9-08	2351683
699 18 733.8-08	1383605	4465277	6,878,342
699 09 770.3-08	1556149	1196757 CH	7,252,803
600 07 636.9-08	1196757	2362590	7,503,997
1297874 CH	1153278	1073505 CH	2024048
1297873 CH			

All specifications are subject to change without notice. Coolheat®, VacRamp™, Dri-Pure®, SampleGuard™, SampleGenie™ and associated logos are trade-marks of Genevac Ltd. Unauthorised use is prohibited.

Your local SP sales representative will be happy to advise on the availability of special order items in addition to the sample holders included in this brochure.

Genevac Ltd.

The Sovereign Centre
Farthing Road
Ipswich
IP1 5AP
United Kingdom

Tel:

Sales: +44 (0) 1473 240000
Service: +44 (0) 1473 243000

Email:

Sales: sales.admin@genevac.co.uk
Service: service@spscientific.com

Application Support: applications@genevac.co.uk

scientificproducts.com

SP Scientific

3538 Main Street
Stone Ridge
NY12484
United States of America

Tel:

Sales: +1 845 267 2211
Service: +1 877 548 4666

Email:

Sales: sales@genevacusa.com
Service: service@spscientific.com

Application Support: applications@genevac.co.uk