


Science with Passion

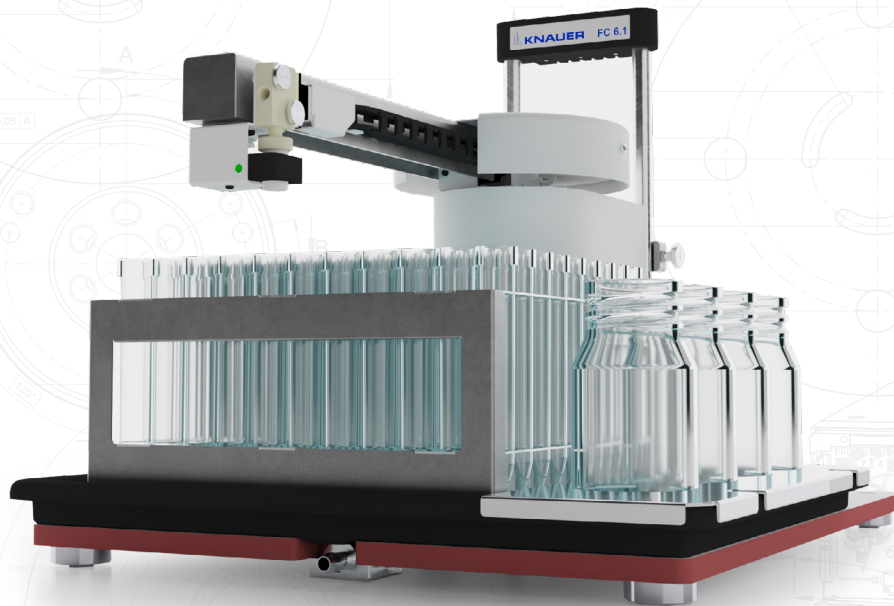

## Fraction Collector FC 6.1

Flexible, reliable and economical

- **Main rack** for tubes and side rack for bottles in one setup
- **Fraction tubes** are highly accessible for sampling
- **Wide flow rate range:** handle drops and flow rates up to 250 ml/min
- **Time-, volume-, drop- or threshold-based** fraction collection
- **Capillary guide** for proper capillary routing
- **Biocompatible** wetted materials
- **Leakage tray** with drain – just in case
- **OEM ready** – communication protocol available

# Fraction collector to handle any collecting challenge

The new FC 6.1 is a small, versatile, and reliable fraction collector. There are two variants available, one designed for aqueous eluents in FPLC and the other for HPLC applications. It can be used with 1/16" or 1/8" tubing for flow rates from 0.1 to 250 ml/min. With its easy-to-change drop formers, the FC 6.1 can be quickly equipped for biocompatibility.

Its main rack and a smaller side rack are accessible from three sides and the top. The adjustable height makes it possible to use bottles of up to 1 l. The FC 6.1 can be placed on the KNAUER AZURA® devices and is supported by PurityChrom® 6, Mobile Control and ClarityChrom software.

## Main Racks for ...

165 tubes, 1/1.5/2 ml each (11 mm)

99 tubes, 15 ml each (17 mm)

30 tubes, 50 ml each (31 mm)

8 round (72 mm) or square (65 mm) bottles, 250 ml each

15 round bottles (56 mm), 100 ml each

## Side Racks for ...

4 round bottle, 100 ml each

3 round/square bottle, 250 ml each

The main racks are not included in the scope of delivery of the FC 6.1 and must be ordered separately.



think **LC**. think **KNAUER**.

Find out more at

<https://www.knauer.net/liquid-handling-fraction-collectors>

(U)HPLC • Prep. LC • FPLC • SMB • LNP • Osmometry

KNAUER Wissenschaftliche Geräte GmbH • Hegauer Weg 38 • 14163 Berlin • +49 30 809727-0 • info@knauer.net • www.knauer.net